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
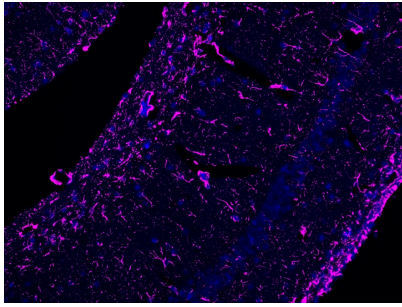
Validation Data

IF Staining using Biotin-TSA Reagent PR30029. FFPE mouse brain tissue was stained with HRP conjugated GFAP monoclonal antibody at 0.5 μ g/mL followed by signal amplification using with Biotin-TSA Reagent PR30029 for 10 minutes at room temperature. Streptavidin-CorLite® Plus 647 Conjugate (Cat.NO. CL647-PF00030) was diluted a... 1:500 and reacted for 10 minutes at room temperature for detection. Cell nucleus (blue) was stained with DAPI.

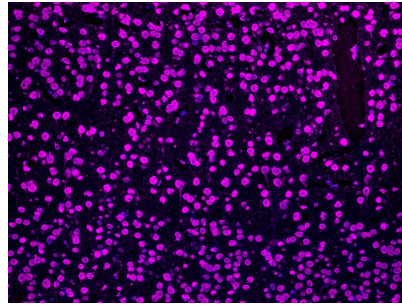IF Staining using Biotin-TSA Reagent PR30029. FFPE mouse brain tissue was stained with HRP conjugated TDP43 rabbit recombinant antibody at 0.5 μ g/mL followed by signal amplification using with Biotin-TSA Reagent PR30029 for 10 minutes at room temperature. Streptavidin-CorLite® Plus 647 Conjugate (Cat.NO. CL647-PF00030) was diluted a... 1:500 and reacted for 10 minutes at room temperature for detection. Cell nucleus (blue) was stained with DAPI.

For technical support and original validation data for this product please contact

T: 027-87531629

E: Proteintech-CN@ptglab.com

W: ptgcn.com

This product is exclusively available under Proteintech Group brand and is not available to purchase from any other manufacturer.